

United States Department of Agriculture  
Center for Veterinary Biologics

Testing Protocol

Fluorescent Antibody Staining Procedure for Detection of Viral Antigens

Date: July 11, 2017  
Number: VIRPRO1014.06  
Supersedes: VIRPRO1014.05, October 10, 2014  
Contact: Alethea M. Fry, (515) 337-7200  
Peg A. Patterson  
Mark A. Troendle

Approvals: /s/Geetha B. Srinivas Date: 20Jul17  
Geetha B. Srinivas, Section Leader  
Virology

United States Department of Agriculture  
Animal and Plant Health Inspection Service  
P. O. Box 844  
Ames, IA 50010

Mention of trademark or proprietary product does not constitute a guarantee or warranty of the product by USDA and does not imply its approval to the exclusion of other products that may be suitable.

|  |
|--|
| Entered into CVB Quality Management System<br>by: <u>/s/Linda S. Snavelly</u> <u>20Jul17</u><br>Linda S. Snavelly Date<br>Quality Management Program Assistant |
|--|

UNCONTROLLED COPY

**Fluorescent Antibody Staining Procedure for Detection of Viral Antigens**

**Table of Contents**

- 1. Introduction**
- 2. Materials**
  - 2.1 Equipment/instrumentation**
  - 2.2 Reagents/supplies**
- 3. Preparation for the Test**
  - 3.1 Personnel qualifications/training**
  - 3.2 Preparation of equipment/instrumentation**
  - 3.3 Preparation of reagents/control procedures**
  - 3.4 Fixation of cell cultures**
- 4. Performance of the Test**
  - 4.1 Direct FA/Indirect FA**
  - 4.2 Reading**
- 5. Interpretation of the Test Results**
- 6. Report of Test Results**
- 7. References**
- 8. Summary of Revisions**

**Appendices**

## Fluorescent Antibody Staining Procedure for Detection of Viral Antigens

### 1. Introduction

This testing protocol (PRO) describes a fluorescent antibody staining method for detection of viral antigens. The protocol uses fluorescein isothiocyanate (FITC) conjugated antibodies for visualization of viral antigen-antibody complexes by ultraviolet (UV) light microscopy. The direct fluorescent antibody (FA) method uses a specific FITC-labeled antibody which binds to the viral antigen directly. The indirect fluorescent antibody (IFA) method uses two specific antibodies: 1) unlabeled primary antibody which binds to a specific viral antigen, and 2) FITC-labeled anti-species secondary antibody which binds to the primary antibody-antigen complex. These methods may be used for:

- Detection of extraneous agents in master seed viruses (MSV)
- Detection of extraneous agent in master cells stocks (MCS)
- Determining MSV neutralization
- Performance of viral titrations
- Performance of identity testing
- Detection of extraneous agents in ingredients of animal origin (FBS. Etc.,)
- Reagent evaluation

### 2. Materials

#### 2.1 Equipment/instrumentation

Equivalent equipment or instrumentation may be substituted for any brand names listed below.

2.1.1 Aerobic incubator,  $36^{\circ}\pm 2^{\circ}\text{C}$ , humidified

2.1.2 Laminar flow biological safety cabinet

2.1.3 Ultraviolet (UV) light microscope

2.1.4 Chemical fume hood

#### 2.2 Reagents/supplies

Equivalent reagents or supplies may be substituted for reagents listed below.

2.2.1 Deionized water (DI)

2.2.2 Acetone (100% or 80% in DI) will be referenced throughout this document; however, other fixatives may be used as appropriate.

**UNCONTROLLED COPY**

**Fluorescent Antibody Staining Procedure for Detection of Viral Antigens**

**2.2.3** Phosphate buffered saline (PBS), 0.01M, pH 7.2, National Centers for Animal Health (NCAH) Media #30054

**2.2.4** FA rinse buffer NCAH Media #30446

**2.2.5** Evan's blue biological stain (EBBS), 0.1% stock solution in DI

**2.2.6** Antigen-specific FITC-conjugated polyclonal antiserum or conjugated monoclonal antibody (direct FA)

**2.2.7** Primary antibody, polyclonal antiserum or monoclonal antibody, specific to viral antigens (IFA)

**2.2.8** FITC-conjugated secondary antibody or anti-species conjugate specific to the species in which the primary antibody was produced (IFA)

**2.2.9** Chambered slides: 1-, 2-, 4-, or 8-well

**2.2.10** Cell culture plates, polystyrene: 24-, 48-, or 96-well

**2.2.11** Microscope slide staining dish with rack or Coplin staining jar

**2.2.12** Miscellaneous lab supplies: beakers, pipettes, wash bottle

**2.2.13** Mounting medium (50% glycerol and 50% FA rinse buffer)

**3. Preparation for the Test**

**3.1 Personnel qualifications/training**

Personnel performing this protocol must have experience or training in aseptic biological laboratory techniques, cell culture, preparation, handling, and disposal of biological agents and chemicals. Personnel must also have knowledge of safe operating procedures and training in the operation of the necessary laboratory equipment listed in **Section 2.1**.

**3.2 Preparation of equipment/instrumentation**

Turn on biological safety cabinet at least 15 minutes before use.

## Fluorescent Antibody Staining Procedure for Detection of Viral Antigens

### 3.3 Preparation of reagents/control procedures

On the day of the FA staining, dilute reagents in **Sections 2.2.6 through 2.2.8** with PBS to a previously determined optimal concentration for the detection of antigens.

### 3.4 Fixation of cell cultures

**Caution: Use fixative in a fume hood or biological safety cabinet. Gloves must be worn to avoid skin contact. Fixatives may be flammable; keep away from sources of heat or flame. Collect used acetone for proper disposal.**

#### 3.4.1 Chamber Slides

1. Decant media from chamber slides into an autoclavable container and remove the plastic walls, leaving slides with the gasket attached. Handle and dispose of all viral fluids according to the current version of **VIRSOP1101**, *Disposal of Live Agents*.

2. Place slides in a slide rack and immerse the loaded slide rack in a staining dish filled with PBS, wash for  $10 \pm 5$  minutes.

**Note: The decision to rinse slides with PBS is optional and is based upon the condition of the cell culture monolayer as observed by the technical staff.**

3. Remove slides from PBS and place in a staining dish filled with acetone for  $10 \pm 5$  minutes. Use 100% acetone for glass slides and 80% acetone for plastic slides. Remove the slides and allow to air dry.

4. If cell preparations are not immediately stained, store monolayers at  $-20^{\circ} \pm 2^{\circ} \text{C}$  or colder until use. Length of acceptable storage will vary for each antigen. Adequacy for use is monitored by FA or IFA staining.

#### 3.4.2 Cell culture plates or dishes

1. Decant the media from plate/dish into an autoclavable container. Handle and dispose of all viral fluids according to the current version of **VIRSOP1101**.

2. With a wash bottle, gently rinse cell monolayer with PBS and decant.

**Note: The decision to rinse slides with PBS is optional and is based upon the condition of the cell culture monolayer as observed by the technical staff.**

**UNCONTROLLED COPY**

### Fluorescent Antibody Staining Procedure for Detection of Viral Antigens

3. If a PBS rinse is used, fill the plate/dish with PBS and allow cell culture layer to soak for  $10 \pm 5$  minutes.
4. Decant PBS from plate/dish.
5. Fix cell monolayer with 80% acetone for  $10 \pm 2$  minutes.
6. Decant acetone into a suitable container; gently blot vessel onto an absorbent surface and allow to air dry.
7. If cell preparations are not immediately stained, store vessels at  $-20^{\circ} \pm 2^{\circ}\text{C}$  or colder until use. Length of acceptable storage will vary for each antigen. Adequacy for use is monitored by FA or IFA staining.

## 4. Performance of the Test

### 4.1 Direct FA/Indirect FA

- 4.1.1 Apply appropriate volume of specific fluorescein-labeled antibody (direct FA) or primary antibody (IFA) to cover the fixed cell monolayer (**Appendix**). Monolayers should not be allowed to dry during staining process.
- 4.1.2 Incubate at  $36^{\circ} \pm 2^{\circ}\text{C}$  for 25 to 65 minutes in a humidified chamber.
- 4.1.3 Decant antibody from monolayers. Rinse at least twice in FA rinse buffer or PBS for  $5 \pm 2$  minutes per rinse. After final rinse, gently tap vessel onto absorbent material to remove excess buffer or PBS buffer.
- 4.1.4 For IFA, repeat **Sections 4.1.1 through 4.1.3** using a secondary FITC-labeled antibody.
- 4.1.5 Gently tap excess FA rinse buffer from monolayers.
- 4.1.6 Monolayers may be rinsed with DI water to remove buffer salts.

**Note: Counterstaining FITC-stained monolayers is often desirable to quench nonspecific background fluorescence or to heighten contrast between specific staining and background. Refer to Section 4.1.7.**

- 4.1.7 (**Optional**) For counterstaining, dilute EBBS stock solution directly in conjugate (1:1000-1200) or in DI (1:1000). Apply EBBS diluted in DI to monolayers and incubate at  $36^{\circ} \pm 2^{\circ}\text{C}$  for  $20 \pm 10$  minutes in a humidified

**UNCONTROLLED COPY**

### Fluorescent Antibody Staining Procedure for Detection of Viral Antigens

chamber. Decant EBBS and quickly rinse with DI. Gently tap vessel onto an absorbent surface.

#### 4.2 Reading

**4.2.1** Monolayers should be read wet and immediately after staining. Vessels are retained for review and shall be stored in the dark at  $4^{\circ} \pm 2^{\circ}\text{C}$  for up to 7 days. Prior to storage, monolayers must be air dried or overlaid with mounting medium. Air dried monolayers must be re-moistened using mounting medium, FA rinse buffer or DI water before review.

**4.2.2** Stained cell cultures are observed at 100X to 200X magnification in a darkened room with the use of a UV light microscope.

**4.2.3** Record results at the time of observation on the appropriate paperwork.

#### 5. Interpretation of the Test Results

**5.1** After staining, each group of monolayers shall be examined for the presence of specific fluorescence attributable to the presence of extraneous viruses. Cells exhibiting specific apple-green fluorescence are considered positive. Non-infected controls should have little nonspecific background fluorescence and be distinguishable from infected controls.

**Note: The above are general guidelines for evaluating stained monolayers. The Agent Contact shall review monolayers to confirm results.**

**5.2** Specifically for MSV testing, a **VALID** test requires:

**5.2.1** The negative control monolayers are free from specific fluorescence. If the negative control monolayers contain specific fluorescence, this indicates possible viral cross contamination, the test is **INVALID**, and must be repeated.

**5.2.2** The positive control monolayers exhibit the expected fluorescence. If specific fluorescence is absent in the positive monolayers, the test is **INVALID**, and must be repeated.

**Note: The positive controls must contain a sufficient number of infected cells to easily determine positive status; however, non-infected cells should also be present in order to differentiate positive cells from negative cells.**

**5.2.3** The MSV inoculated cells are free from specific fluorescence to the MSV. If the MS inoculated cells contain specific fluorescence to the MSV, this indicates

**UNCONTROLLED COPY**

### Fluorescent Antibody Staining Procedure for Detection of Viral Antigens

incomplete neutralization and/or replication of MSV, the test is INVALID, and must be repeated.

**5.2.4** If the test is valid and the MSV test monolayers are equivocal, then the test is declared INCONCLUSIVE, and must be repeated.

**5.2.5** If the test is valid and the MSV test monolayers are negative for specific fluorescence, then the MSV is SATISFACTORY.

**5.2.6** If the test is valid and specific fluorescence is found in the MSV test monolayers, then the test MSV is UNSATISFACTORY. At supervisory discretion, the test may be repeated to confirm result.

## 6. Report of Test Results

All records are kept in accordance with the current recordkeeping practices. Test results will be reviewed and signed by knowledgeable staff (such as the Agent Contact, Microbiologist, or Lead Technician). Test results are then entered into the current reporting system and released to the reviewer for distribution to the firm.

## 7. References

Title 9, *Code of Federal Regulations*, sections 113.47 and 113.309, U.S. Government Printing Office, Washington, DC.

## 8. Summary of Revisions

### Version .06

- The FA use list has been updated to include reagents used to detect and test for Seneca Valley virus.

### Version .05

- The Contact information has been updated; however, the Virology Section has elected to keep the same next review date for the document.

### Version .04

- The Contact information has been updated.

**UNCONTROLLED COPY**



**Fluorescent Antibody Staining Procedure for Detection of Viral Antigens**

- **1:** The FA use list has been expanded to include MCS and ingredients of animal origin.
- **2.2:** Redundant equipment has been removed from the list.
- **3.4:** An optional PBS wash has been included.
- **4.1.2:** A time range for the incubation time of fluorescein-labeled antibody or primary antibody has been added.
- **4.1.3:** The antibody rinse of slides or plates has been changed from 3 to 2. This reflects the common practice which is in place at the CVB. Also added the option of using PBS to rinse plates or slides.
- **4.1.6:** The option of using a DI rinse to remove residual salts has been added.
- **4.2.1:** The option of using wetting slides or plates with either PBS or DI water, prior to reading has been added.
- **6:** The description of staff able to review results has been updated.
- **Appendix I:** Volume ranges are now included.
- **Appendix II:** A table which includes the recommended positive control viruses, and their use dilutions, and the test detection reagents and applicable use dilutions has been added to the document.

**Version .03**

- The Contact information has been updated.
- National Veterinary Services Laboratories (NVSL) has been changed to National Centers for Animal Health (NCAH) throughout the document.

**Version .02**

- The title of this document has been revised and the document has been rewritten to superseded VIRSOP1500 and VIRPRO1010.
- An appendix has been added to include volumes of reagents used for various vessels.

**UNCONTROLLED COPY**

Fluorescent Antibody Staining Procedure for Detection of Viral Antigens

**Appendix I**  
**Volume of Antibody to be used for FA Staining per Vessel**

| <b>Vessel</b>    | <b>Volume/chamber or well (<math>\mu</math>L)</b> |
|------------------|---|
| 2-chamber slides | 200-300   |
| 4-chamber slides | 100-150   |
| 8-chamber slides | 50-75   |
| 24-well plates   | 200-300   |
| 48-well plates   | 100-150   |
| 96-well plates   | 50-75   |
| 60 mm plates     | 2000  |

Fluorescent Antibody Staining Procedure for Detection of Viral Antigens

Appendix II

| Agent                                   | Detection Method   | Primary antibody | Concentration | Conjugate identification  | Conjugate concentration | Cell line         | Required by 9CFR | Positive control (PC)        | Lot Titer              | PC concentration (log) |
|---|--------------------|------------------|---------------|---------------------------|-------------------------|-------------------|------------------|------------------------------|------------------------|------------------------|
| BVDV-1                                  | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-BVD-10ML             | <i>Undiluted</i>        | BT                | √                | 00-17                        | 10 <sup>6.6</sup> /2mL | 3-4 (1:1000)           |
| BVDV-2                                  | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-BVD-10ML             | <i>Undiluted</i>        | BT                | √                | 05-15                        | 10 <sup>6.3</sup> /2mL | 3-4 (1:1000)           |
| Blue Tongue Serotype 10                 | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-FA-BTV-MAB-10ML        | <i>Undiluted</i>        | BT                | √                | 85-9                         | 10 <sup>4.5</sup> /1mL | 2.5 (1:100)            |
| Rabies                                  | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | <i>Fujireloio 800-092</i> | <i>1:100</i>            | <i>Any</i>        | √                | <i>Virology Slides</i>       | <i>N/a</i>             | <i>N/a</i>             |
| Reovirus                                | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-REO-10ML             | <i>Undiluted</i>        | Vero              | √                | 01-17                        | 10 <sup>5.0</sup> /1mL | 3 (1:100)              |
| <b>Bovine</b>                           |                    |                  |               |                           |                         |                   |                  |                              |                        |                        |
| Bovine adenovirus type 1                | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-BAV1-10ML            | <i>Undiluted</i>        | MDBK-A            | √                | 06-39                        | 10 <sup>5.7</sup> /2mL | 3 (1:100)              |
| Bovine adenovirus type 5                | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-BAV5-10ML            | <i>Undiluted</i>        | BT                | √                | 04-0401                      |                        | 2-3 (1:100)            |
| Bovine coronavirus                      | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-BCV-10ML             | <i>Undiluted</i>        | MDBK              | N/a              | 05-01                        |                        | 2                      |
| Bovine parvovirus                       | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-BPV-10ML             | <i>Undiluted</i>        | BT                | √                | 99-09                        | 10 <sup>5.2</sup> /1mL | 4 (1:10)               |
| Bovine respiratory syncytial virus      | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-BRSV-10ML            | <i>Undiluted</i>        | BT                | √                | 88-16                        | 10 <sup>5.4</sup> /2mL | 3 (1:100)              |
| Infectious bovine rhinotracheitis virus | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-IBR-10ML             | <i>Undiluted</i>        | BT                | N/a              | 84-5, 92-12                  | 10 <sup>6.2</sup> /1mL | 4 (1:100)              |
| Parainfluenza virus type 3              | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-PI#-10ML             | <i>Undiluted</i>        | BT                | N/a              | 84-6                         | 10 <sup>7.0</sup> /1mL | 4 (1:1000)             |
| <b>Porcine</b>                          |                    |                  |               |                           |                         |                   |                  |                              |                        |                        |
| Swine adenovirus                        | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-PAV-10ML             | <i>Undiluted</i>        | <i>B-ST/PK-15</i> | √                | <i>Strain NADC lot 85-17</i> | 10 <sup>6.3</sup> /1mL | 3-4 (1:100-1:1000)     |
| Porcine parvovirus                      | <i>FA (Direct)</i> | <i>N/a</i>       | <i>N/a</i>    | CJ-F-PPV-10ML             | <i>Undiluted</i>        | <i>B-ST/PK-15</i> | √                | <i>Strain NADL lot 87-6</i>  | 10 <sup>4.5</sup> /1mL | 2-3 (1:100)            |

UNCONTROLLED COPY

Fluorescent Antibody Staining Procedure for Detection of Viral Antigens

| Agent  | Detection Method | Primary antibody | Concentration | Conjugate identification               | Conjugate concentration                   | Cell line  | Required by 9CFR | Positive control (PC) | Lot Titer              | PC concentration (log) |
|--|------------------|------------------|---------------|--|---|------------|------------------|-----------------------|------------------------|------------------------|
| Transmissible gastroenteritis virus              | FA (Direct)      | N/a              | N/a           | CJ-F-TGE-10ML                          | Undiluted                                 | B-ST/PK-15 | √                | Strain M59S, lot M59S |                        | 1.5                    |
| Porcine hemagglutinating encephalomyelitis virus | FA (Direct)      | N/a              | N/a           | CJ-F-PHEV-10 ML                        | Undiluted                                 | B-ST/PK-15 | √                | HEV e lot 01-16       |                        | 2                      |
| Porcine circovirus type 1                        |                  |                  |               |  |   |            | N/a              |                       |                        |                        |
| Porcine circovirus type 2                        | FA (Direct)      | N/a              | N/a           | CJ-F-PCV2-10ML                         | Undiluted                                 | B-ST/PK-15 | N/a              | PCV2/6 88 12-4-01     |                        | 2.5-3                  |
| Porcine respiratory and reproductive virus       |                  |                  |               |  |   | Marc 104   | N/a              |                       |                        |                        |
| Swine influenza Virus                            |                  |                  |               |  |   | MDCK       | N/a              |                       |                        |                        |
| Seneca Valley Virus                              | FA (Indirect)    | 88-3669S 4A      | 1:20          | KPL Goat anti-swine Catalog # 02-14-06 | 1:100                                     | CEF        | N/a              | SVV150 31348          | TBD                    | TBD                    |
| <b>Equine</b>                                    |                  |                  |               |  |   |            |                  |                       |                        |                        |
| Equine herpesvirus type 1                        | FA (Direct)      | N/a              | N/a           | CJ-F-ERV-10ML                          | Undiluted                                 | EQ Der     | √                | 5-85                  |                        | 4 (1:10)               |
| Equine herpesvirus type 4                        | FA (Direct)      | N/a              | N/a           | CJ-F-ERV-10ML                          | Undiluted                                 | EQ Der     | √                | 01-03                 | 10 <sup>4.8</sup> /2mL | 3 (1:100)              |
| Equine herpesvirus type 4 (confirmation)         | FA (Indirect)    |                  |               |  |   | EQ Der     |                  |                       |                        |                        |
| Equine arteritis virus                           | FA (Indirect)    | CVB              | 1:100         | 92-10                                  | Secondary: KPL anti-equine (1:100)        | EQ Der     | √                | 94-10                 | 10 <sup>6.1</sup> /2mL | 5 (1:10)               |
| Equine infectious anemia virus                   |                  |                  |               |  |   | EQ Der     | N/a              |                       |                        |                        |
| Equine influenza virus                           | FA (Indirect)    | CVB              | 1:50-1:100    | 06-31                                  | Secondary: NVSL anti-chicken 291 FA 860 2 | MDCK       | N/a              | 92-7 or 1-86 or 79    | Vary                   | 1:100                  |

UNCONTROLLED COPY

Fluorescent Antibody Staining Procedure for Detection of Viral Antigens

| Canine                                   |                      |                    |                    |   |  |             |     |                      |                         |                             |
|--|----------------------|--------------------|--------------------|---|--|-------------|-----|----------------------|-------------------------|-----------------------------|
| Canine coronavirus                       | <i>FA (Direct)</i>   | <i>N/a</i>         | <i>N/a</i>         | CJ-F-CCV-10ML                               | <i>Undiluted</i>                       | <i>CrFK</i> | √   | 07-06<br>or<br>91-28 | 10 <sup>5.2</sup> /1mL  | 3 (1:100)                   |
| Canine distemper virus                   | <i>FA (Direct)</i>   | <i>N/a</i>         | <i>N/a</i>         | CJ-F-CDV-10Ml or<br>CJ-F-CDV-MAB-<br>10ML   | <i>Undiluted</i>                       | <i>MDCK</i> | √   | 99-05                | 10 <sup>4.7</sup> /1mL  | 2.7<br>(1:100)              |
| Canine parvovirus                        | <i>FA (Direct)</i>   | <i>N/a</i>         | <i>N/a</i>         | CJ-F-CPV-MAB-<br>10Ml or CJ-F1-CPV-<br>10ML | <i>Undiluted</i>                       | <i>CrFK</i> | √   | 96-05                | 10 <sup>6.4</sup> /1mL  | 4.4<br>(1:100)              |
| Canine parainfluenza virus               | <i>FA (Direct)</i>   | <i>N/a</i>         | <i>N/a</i>         | CJ-F-CPI-10ML                               | <i>Undiluted</i>                       | <i>Vero</i> | N/a | 96-16                | 10 <sup>8.5</sup> /1mL  | 3-4<br>(1:1000-<br>1:10000) |
| Canine herpes virus                      | <i>FA (Direct)</i>   | <i>N/a</i>         | <i>N/a</i>         | CJ-F-CHV-10ML                               | <i>Undiluted</i>                       | <i>MDCK</i> | N/a | VR552<br>or 97-16    | 10 <sup>6.79</sup> /1mL | 5 (1:100)                   |
| Canine adenovirus type 2                 | <i>FA (Direct)</i>   | <i>FA (Direct)</i> | <i>FA (Direct)</i> | CJ-F-CAV-10mL                               | <i>Undiluted</i>                       | <i>MDCK</i> | N/a | 06-44                | 10 <sup>7.74</sup> /1mL | 5 (1:1000)                  |
| Infectious canine hepatitis virus (CAV1) | <i>FA (Direct)</i>   | <i>FA (Direct)</i> | <i>FA (Direct)</i> | CJ-F-CAV-10ML                               | <i>Undiluted</i>                       | <i>MDCK</i> | N/a | 06-44                | 10 <sup>7.74</sup> /1mL | 5<br>(1:1000)               |
| Canine influenza virus                   |                      |                    |                    |   |  | <i>MDCK</i> | N/a |                      |                         |                             |
| Feline                                   |                      |                    |                    |   |  |             |     |                      |                         |                             |
| Feline infectious peritonitis virus      | <i>FA (Direct)</i>   | <i>N/a</i>         | <i>N/a</i>         | CJ-F-FIP-10ML                               | <i>Undiluted</i>                       | <i>CrFK</i> | √   | 06-42                |                         | 4                           |
| Feline parvovirus                        | <i>FA (Direct)</i>   | <i>N/a</i>         | <i>N/a</i>         | CJ-F-FPL-10ML                               | <i>Undiluted</i>                       | <i>CrFK</i> | √   | 02-14                | 10 <sup>6.83</sup> /1mL | 3<br>(1:1000)               |
| Feline calicivirus                       | <i>FA (Indirect)</i> | <i>CVB</i>         | <i>1:80</i>        | <i>96-02</i>                                | <i>Secondary: KPL anti-cat (1:100)</i> | <i>CrFK</i> | N/a | 96-14                | 10 <sup>5.2</sup> /1mL  | 3 (1:100)                   |
| Feline rhinotracheitis virus             | <i>FA (Direct)</i>   | <i>N/a</i>         | <i>N/a</i>         | <i>CJ-F-FVR-10ML</i>                        | <i>Undiluted</i>                       | <i>CrFK</i> | N/a | 96-15                | 10 <sup>8.26</sup> /1mL | 5<br>(1:1000)               |
| Feline leukemia virus                    | <i>FA (Indirect)</i> | 210-46-<br>FeLV1   | <i>Undiluted</i>   | AB1-FELV and AB2-<br>FELV                   | <i>Undiluted</i>                       | <i>CrFK</i> | N/a | 95-09<br>(Challenge) | 10 <sup>6.0</sup> /1mL  | 4 (1:100)                   |

UNCONTROLLED COPY